



Lonza

Triangle Research Labs – Now Part of Lonza

Triangle Research Labs, LLC | www.TRLcells.com

U.S. Scientific Support: 800 521 0390 | scientific.support@lonza.com

EU/ROW Scientific Support: +32 87 321 611 | scientific.support.eu@lonza.com

Completing Media for Use with Primary Hepatocytes Protocol

- 1) To complete the hepatocyte plating media (MP100):**
 - a. Pour the entire contents of the vial labeled 'Plating Supplement' into the media.
 - b. Sterile filter the completed media

- 2) To complete the hepatocyte maintenance media (MM250):**
 - a. Pour the entire contents of the vial labeled 'Maintenance Supplement' into the media.
 - b. Sterile filter the completed media